

# THE BAKER COMPANY

161 Gatehouse Rd., Sanford, Maine 04073 USA  
"Creating Immaculate Atmospheres"

## SG3 AUX WIRING OPTIONS

**In response to customer requests and effective 08/30/05, new wiring configurations will now be available on the SterilChemGARD's & SterilGARD's. The 0.1 Amps "C" Contact & the Auxiliary Junction Box Options will no longer be available.**

### **SterilChemGARD SG403TX/603TX**

All SG403TX/603TX units will have the Blower Switch Status Wiring Package (333E548) installed as a standard configuration. This package includes wiring and a junction box that connects to the Blower Switch Status relay. This wiring allows the customer to connect the unit to a system to monitor the switch status or to connect to a control circuit for an external blower or damper in their exhaust system.

When the customer wants to monitor additional conditions in the unit, they can order the optional Cabinet Monitor Wiring Package (333E549). This option will include the wiring and hardware for the Blower Switch Status Wiring Package (333E548) plus wiring for connection to the Blower Operational Status relay and the Sash Alarm Status relay. They can choose to connect to one or more of these relays for monitoring or control of external systems.

To upgrade older units, the customer would need to purchase a replacement Membrane Switch Control Board (333E550) that provides the status relays and either the Blower Switch Status Wiring Package (333E548) or the Cabinet Monitor Wiring Package (333E549).

### **SterilGARD**

There will be two options available for the SG403A & SG603A

1. The Blower Switch Status Wiring Package (333E548) will provide wiring and a junction box with connections for the Blower Switch Status relay. The customer will be able to use this wiring to connect to a system to monitor the switch status or to connect to a control circuit for an external blower or damper in the exhaust system.
2. The Cabinet Monitor Wiring Package (333E549) will provide wiring and a junction box with connections for the Blower Switch Status relay, Blower Operational Status relay and Sash Alarm Status relay. The customer can choose to connect to one or more of these relays for monitoring or control of external systems.

161 Gatehouse Rd., Sanford, Maine 04073 USA · (207) 324-8773 · 1-800-992-2537 · FAX (207) 324-2632  
Internet address: [http:// www.bakerco.com](http://www.bakerco.com) or e-mail: [bakerco@bakerco.com](mailto:bakerco@bakerco.com)

*"Creating Immaculate Atmospheres"*